


Laboratory Location (Building/Room):

Completed By:


Date:

**BSC Workspace:**

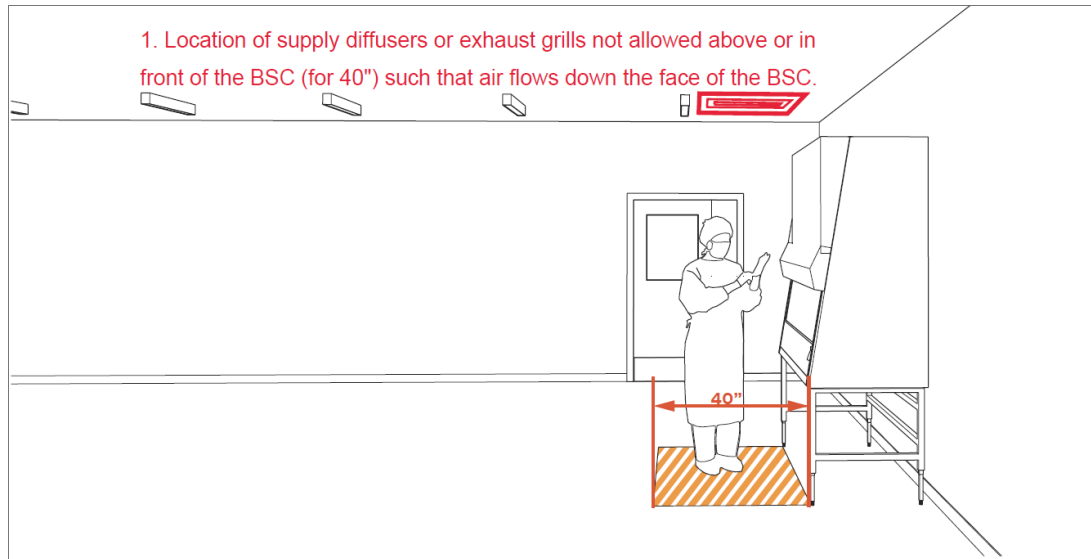

2. Maintain an undisturbed space of 40" around the BSC. \*Undisturbed means free of both permanent (walls, cabinets, equipment) and temporary (door swings, people traffic, cabinet drawers/doors) encroachment.

**Distance to Adjacent Walls & Columns:**

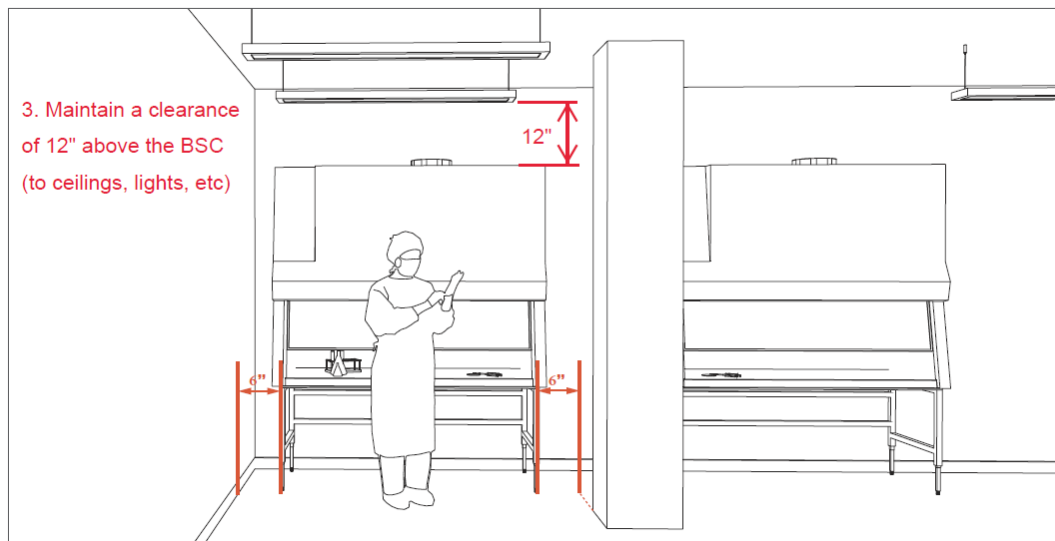

4. Maintain a clearance of 6" on sides and behind the BSC (to walls, columns, benchtops, equipment, etc)


I confirm that this replacement has the appropriate clearances.

I confirm that this replacement does not have the appropriate clearances. I will discuss with DOHS.

I confirm that this replacement has the appropriate clearances.


I confirm that this replacement does not have the appropriate clearances. I will discuss with DOHS.

### BSC Placement Along Opposing Walls:

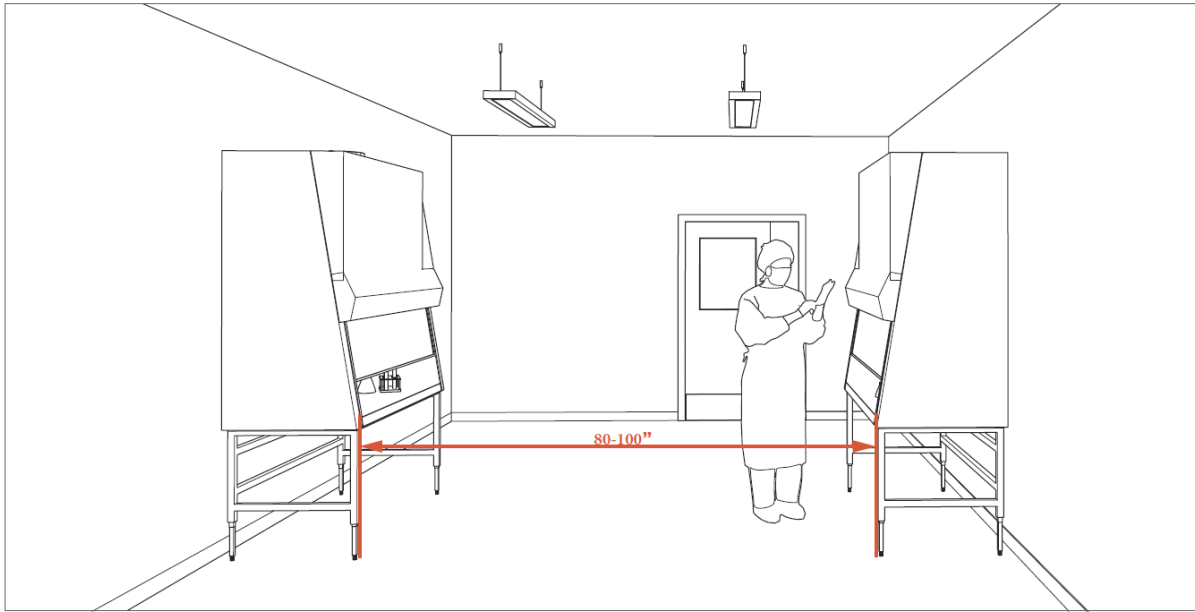

Maintain a distance of 80" between BSCs on opposing walls where there is no potential traffic between BSCs and 100" when there is potential traffic.

### Distance to Opposing Bench Tops:

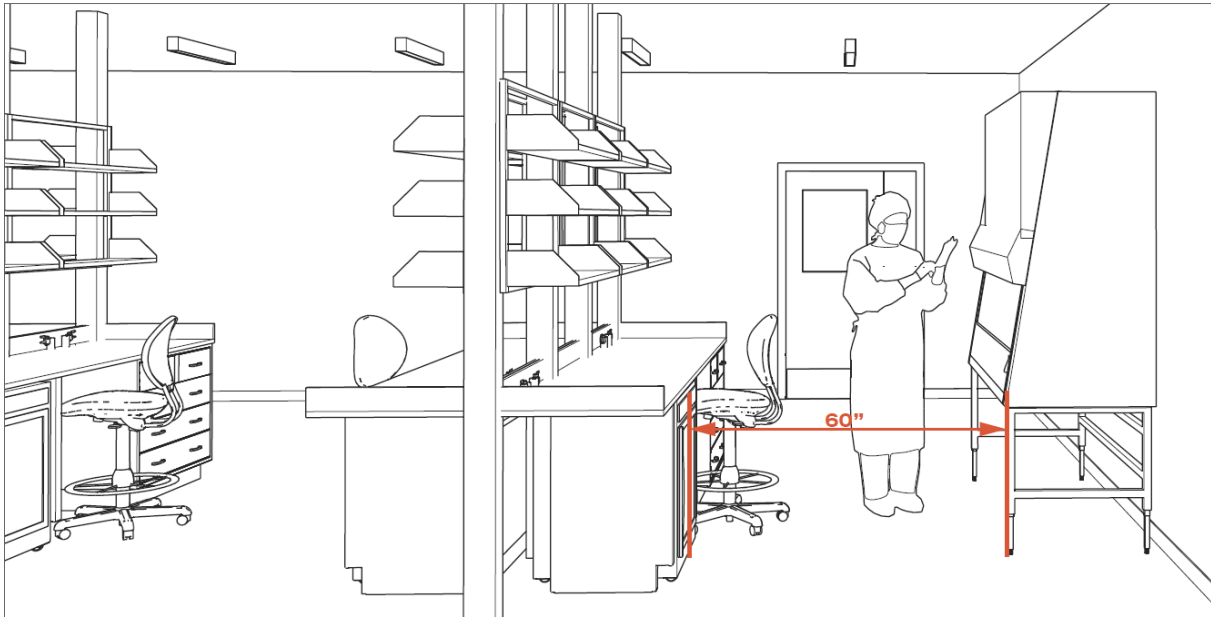

Maintain a clearance of 60" in front of the BSC if there is any possibility of traffic behind the occupant working at the BSC.

I confirm that this replacement has the appropriate clearances.

I confirm that this replacement does not have the appropriate clearances. I will discuss with DOHS.

Not Applicable

I confirm that this replacement has the appropriate clearances.

I confirm that this replacement does not have the appropriate clearances. I will discuss with DOHS.

Not Applicable